


# Viewcen

## Introduction

The **ViewCen** utility provides a quick method for visually checking the quality of the raw camera data stored in a **.cen** file. First select the files you wish to view in the **Input files** list box of the AMASS shell, and then click on the **ViewCen** function button.

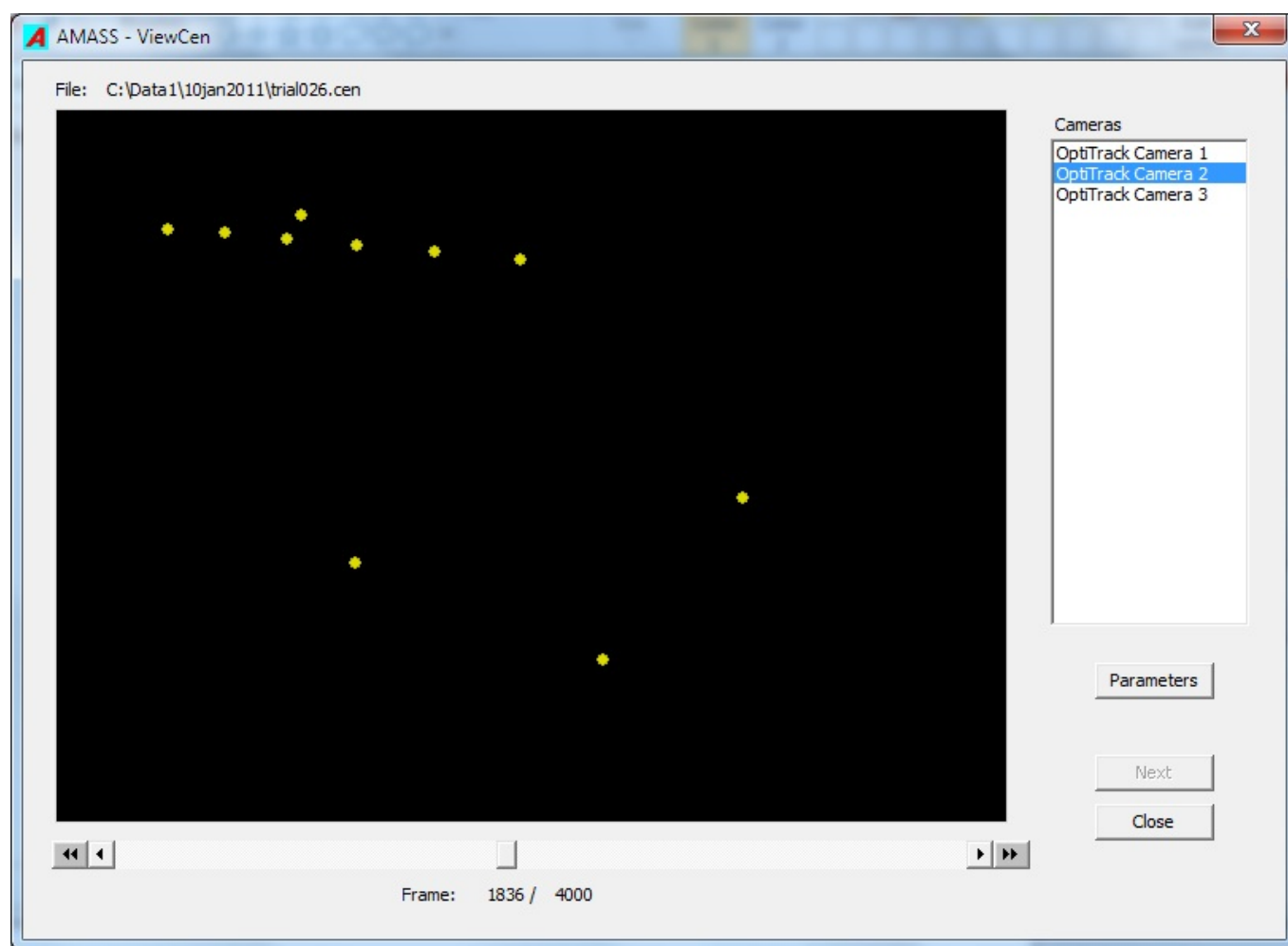

The image area shows the full view as seen by the selected camera, and the slider may be used to access any frame. Clicking a double arrow button causes the frames to be replayed at the collected frame rate. Clicking the button a second time doubles the replay rate, etc. Click on the slider area to stop the replay. The keyboard arrow keys may also be used to change the camera selection.


**ViewCen** will run automatically after each raw camera file conversion if the **ViewCen** box was checked in the **Input** page of the Project parameters.

**ViewCen** may also be run as a standalone program.

## Button functions

### Parameters

This button brings up a window displaying some internal parameters for the `.cen` file.

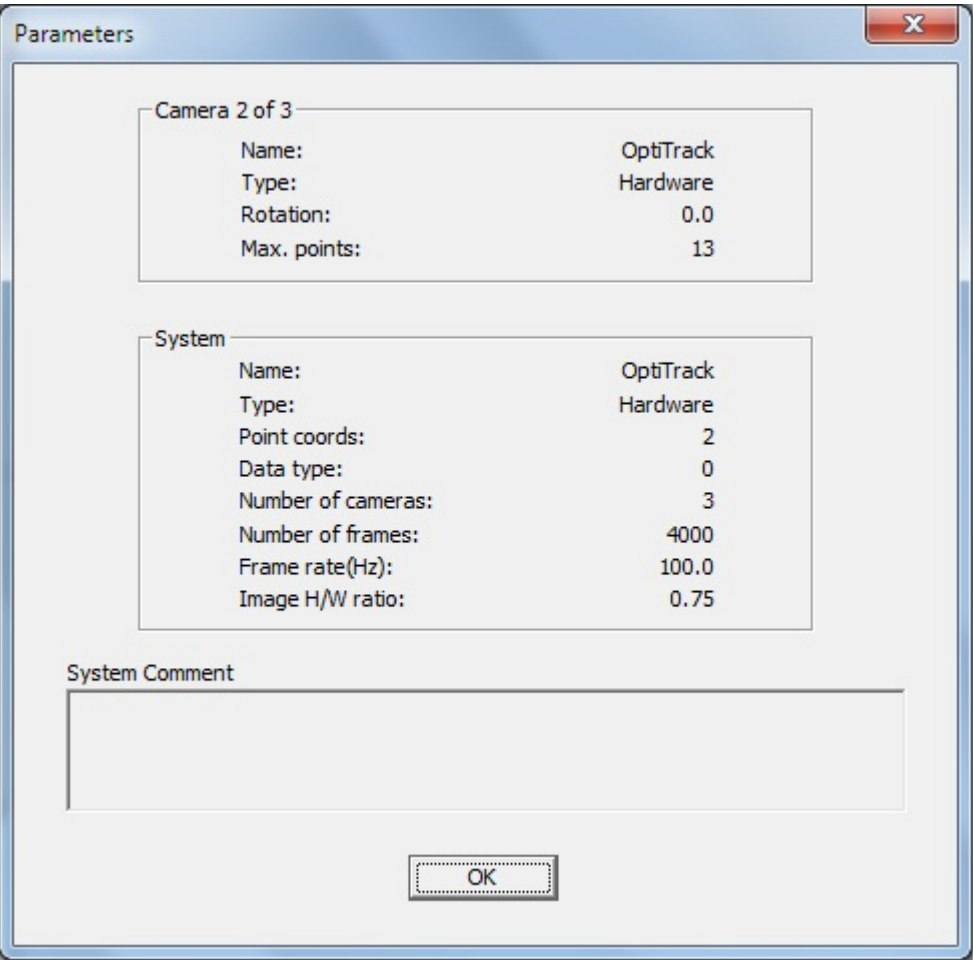


### Next

Displays the next `.cen` file if more than one `.cen` file was selected before the **ViewCen** button was clicked in the AMASS shell.

### Close

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Use this button to exit the **ViewCen** program.



From:

<https://has-motion.com/wiki/> - **Wiki Documentation Site**

Permanent link:

<https://has-motion.com/wiki/doku.php?id=other:amass:documentation:viewcen&rev=1721230911>

Last update: **2024/07/17 15:41**

