

.cen files]] will be recorded can be set by clicking 'browse...'. capture_browse.jpg if capture is launched by amass, capture will use the [[other:report3d:documentation:centroid_files|.cen file]] path the user has set in amass. if capture is launch by the user manually (i.e., double-clicking the capture icon on the desktop), then capture will use: - the last directory set by the user within capture or - the application folder if a directory has never been selected by the user. a separate edit box contains the name of the file. after each .cen file is captured and saved to disk the program will automatically append (or increment) a number at the end of the filename. for example, trail.cen will changed to trail2.cen, etc. if the [[other:report3d:documentation:centroid_files|.cen extension]] is omitted by the user, it will be added automatically to the filename. } } }

From:

<https://has-motion.com/wiki/> - **Wiki Documentation Site**

Permanent link:

https://has-motion.com/wiki/doku.php?id=other:capture2d:connecting_to_cameras&rev=1718804089

Last update: **2024/06/19 13:34**

