


Playback

Once the trial is complete, Capture enters playback mode, so that the captured images can be reviewed.

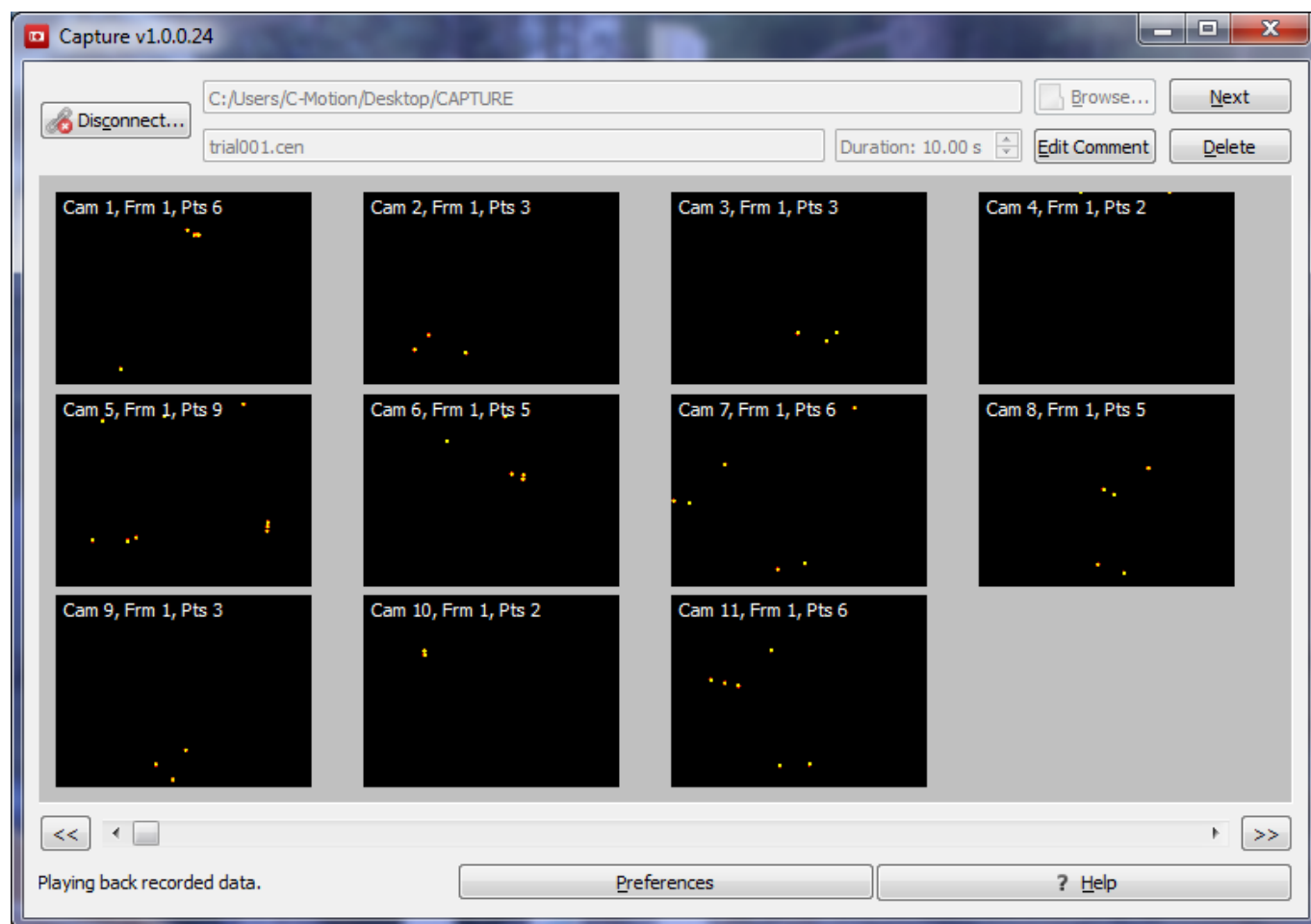

The [.cen files](#) can be viewed by dragging the scroll bar or by clicking the double arrow button. Comments can be added to [.cen files](#) with the 'Edit Comment' button before or after the recording. The trial will be saved to the chosen directory if the 'Next' button is clicked, or the file can be deleted. Capture will then return to the preview mode, and data can be collected again. If the previous trial was saved, the filename will auto-increment to prevent data from being overwritten.

From:
<https://has-motion.com/wiki/> - Wiki Documentation Site

Permanent link:
<https://has-motion.com/wiki/doku.php?id=other:capture2d:playback&rev=1721228966>

Last update: 2024/07/17 15:09

