

once the start recording button is pressed a countdown will begin as indicated by the status bar at the bottom of the dialog and then the data collection will start. when recording starts the status bar then indicates the progress of the recording and the number of frames recorded. capture_trails.png in this figure the marker trails have been turned on. this is a useful feature when capturing the calibration trial because it provides a quick overview of the coverage of the wand through the view of each camera. in this example camera_12 will likely not calibrate well because the wand has not covered a sufficient region of the viewer

From:

<https://has-motion.com/wiki/> - **Wiki Documentation Site**

Permanent link:

https://has-motion.com/wiki/doku.php?id=other:capture2d:recording_data&rev=1718801144

Last update: **2024/06/19 12:45**

