


DSX_Configuration

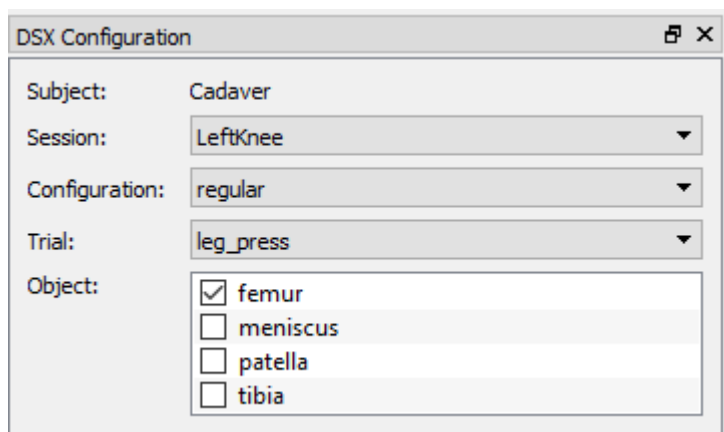

DSX Configuration

Subject: Cadaver

Session: LeftKnee

Configuration: regular

Trial: leg_press

Object:

- ☒ femur
- ☐ meniscus
- ☐ patella
- ☐ tibia

- **Subject** This field is automatically populated with the ID of the subject when you load a subject file.
- **Session** This drop down menu is automatically populated with all of the sessions available for the current subject.
- **Configuration** This drop down menu is automatically populated with all of the configurations available for the current session.
- **Trial** This drop down menu is automatically populated with all of the trials available for the current configuration.
- **Object** This list is automatically populated with all of the objects in the current trial. When an object is selected, its POIs are displayed in the *Points of Interest* widget.

From:
<https://has-motion.com/wiki/> - Wiki Documentation Site

Permanent link:
https://has-motion.com/wiki/doku.php?id=other:dsx:locate3d:dsx_configuration&rev=1721228970

Last update: 2024/07/17 15:09

