


DSX Configuration

X4D's DSX Configuration widget allows users to view and modify settings relevant to the whole DSX Suite, namely information related to subjects, sessions, and trial.

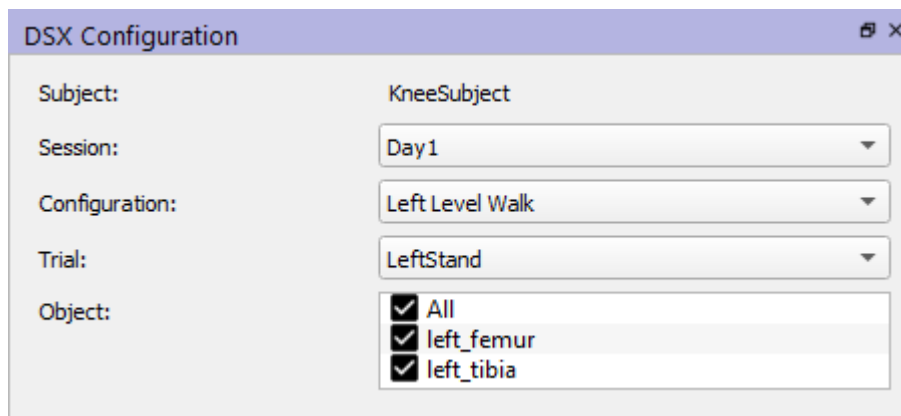

The screenshot shows a window titled "DSX Configuration" with a close button in the top right corner. Inside the window, there are five labeled fields on the left and their corresponding values on the right:

- Subject:** KneeSubject
- Session:** Day1 (dropdown menu)
- Configuration:** Left Level Walk (dropdown menu)
- Trial:** LeftStand (dropdown menu)
- Object:** A list with three items, each preceded by a checked checkbox:
 - All
 - left_femur
 - left_tibia

- **Subject** This field is automatically populated with the ID of the subject when you load a subject file.
- **Session** This drop down menu is automatically populated with all of the sessions available for the current subject.
- **Configuration** This drop down menu is automatically populated with all of the configurations available for the currently selected session.
- **Trial** This drop down menu is automatically populated with all of the trials available for the currently selected configuration.
- **Object** This list is automatically populated with all of the objects in the currently selected trial. Each object is enabled if it is trackable, meaning that it has a 3D image file associated with it, for generating DRRs.

From:
<https://has-motion.com/wiki/> - **Wiki Documentation Site**

Permanent link:
https://has-motion.com/wiki/doku.php?id=other:dsx:x4d:dsx_configuration

Last update: **2025/06/02 17:53**

