


DSX Configuration

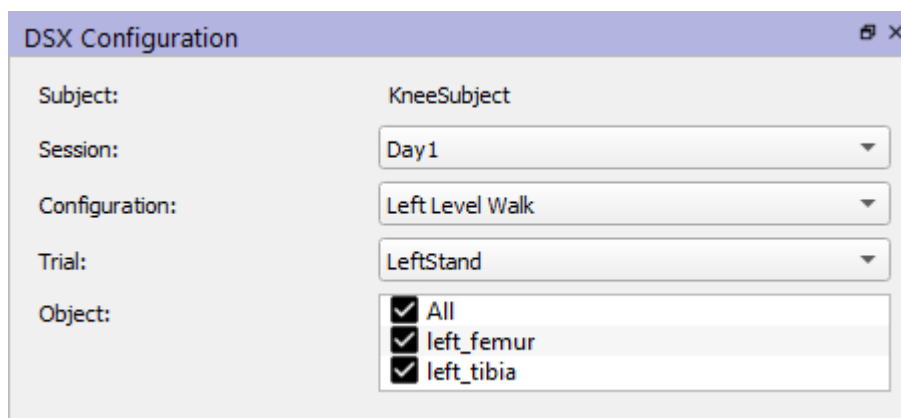

Subject:	KneeSubject
Session:	Day1
Configuration:	Left Level Walk
Trial:	LeftStand
Object:	☒ All ☒ left_femur ☒ left_tibia

- **Subject** This field is automatically populated with the ID of the subject when you load a subject file.
- **Session** This drop down menu is automatically populated with all of the sessions available for the current subject.
- **Configuration** This drop down menu is automatically populated with all of the configurations available for the currently selected session.
- **Trial** This drop down menu is automatically populated with all of the trials available for the currently selected configuration.
- **Object** This list is automatically populated with all of the objects in the currently selected trial. Each object is enabled if it is trackable, meaning that it has a 3D image file associated with it, for generating DRRs.

From:
<https://has-motion.com/wiki/> - **Wiki Documentation Site**

Permanent link:
https://has-motion.com/wiki/doku.php?id=other:dsx:x4d:dsx_configuration&rev=1721230983

Last update: **2024/07/17 15:43**

