

IORgait - Marker Placement

The IOR marker set is comprised of 36 markers. It is best if marker placement done by an individual who is experienced with palpating anatomical positions.

The anatomical location of each target is referenced to the description in the [Color Atlas of Skeletal Landmark Definitions](#). This text is useful for further description of target location as well as recommendations for palpation.

[illegible]

IORfoot - Marker Placement

The anatomical location of each target is referenced to the description in the [Color Atlas of Skeletal Landmark Definitions](#). This text is useful for further description of target location as well as recommendations for palpation.

The FMB and SMB target locations are not described in the [Color Atlas](#), but are described in the [Leardini multi-segment foot paper](#).

Marker Measurement Listed by from knees to feet	Marker Name	Description	Reference Page	URFXF1	Proximal tip of the lateral head of the tibia	URFXF2	Most anterior border of the tuberosity	URFXF3	Lateral border of the lateral malleolus	URFXF4	Most medial border of the lateral malleolus	URFXF5	Aspect of the lateral border of the calcaneus	URFXF6	Sustentaculum of the Calcaneus	URFXF7	Lateral border of the second metatarsal	URFXF8	Navicular tuberosity	URFXF9	Tuberosity of 5th Metatarsal	URFXF10	Base of First Metatarsal	URFXF11	Base of Second Metatarsal	URFXF12	Dorsal margin of the first metatarsal	URFXF13	Dorsal aspect of the second metatarsal	URFXF14	Dorsal margin of the third metatarsal	URFXF15	Dorsal margin of the fourth metatarsal	URFXF16	Dorsal margin of the fifth metatarsal	URFXF17	Dorsal margin of the first phalanx	URFXF18	Dorsal margin of the second phalanx	URFXF19	Dorsal margin of the third phalanx	URFXF20	Dorsal margin of the fourth phalanx	URFXF21	Dorsal margin of the fifth phalanx	URFXF22	Dorsal margin of the first phalanx	URFXF23	Dorsal margin of the second phalanx	URFXF24	Dorsal margin of the third phalanx	URFXF25	Dorsal margin of the fourth phalanx	URFXF26	Dorsal margin of the fifth phalanx	URFXF27	Dorsal margin of the first phalanx	URFXF28	Dorsal margin of the second phalanx	URFXF29	Dorsal margin of the third phalanx	URFXF30	Dorsal margin of the fourth phalanx	URFXF31	Dorsal margin of the fifth phalanx	URFXF32	Dorsal margin of the first phalanx	URFXF33	Dorsal margin of the second phalanx	URFXF34	Dorsal margin of the third phalanx	URFXF35	Dorsal margin of the fourth phalanx	URFXF36	Dorsal margin of the fifth phalanx	URFXF37	Dorsal margin of the first phalanx	URFXF38	Dorsal margin of the second phalanx	URFXF39	Dorsal margin of the third phalanx	URFXF40	Dorsal margin of the fourth phalanx	URFXF41	Dorsal margin of the fifth phalanx	URFXF42	Dorsal margin of the first phalanx	URFXF43	Dorsal margin of the second phalanx	URFXF44	Dorsal margin of the third phalanx	URFXF45	Dorsal margin of the fourth phalanx	URFXF46	Dorsal margin of the fifth phalanx	URFXF47	Dorsal margin of the first phalanx	URFXF48	Dorsal margin of the second phalanx	URFXF49	Dorsal margin of the third phalanx	URFXF50	Dorsal margin of the fourth phalanx	URFXF51	Dorsal margin of the fifth phalanx	URFXF52	Dorsal margin of the first phalanx	URFXF53	Dorsal margin of the second phalanx	URFXF54	Dorsal margin of the third phalanx	URFXF55	Dorsal margin of the fourth phalanx	URFXF56	Dorsal margin of the fifth phalanx	URFXF57	Dorsal margin of the first phalanx	URFXF58	Dorsal margin of the second phalanx	URFXF59	Dorsal margin of the third phalanx	URFXF60	Dorsal margin of the fourth phalanx	URFXF61	Dorsal margin of the fifth phalanx	URFXF62	Dorsal margin of the first phalanx	URFXF63	Dorsal margin of the second phalanx	URFXF64	Dorsal margin of the third phalanx	URFXF65	Dorsal margin of the fourth phalanx	URFXF66	Dorsal margin of the fifth phalanx	URFXF67	Dorsal margin of the first phalanx	URFXF68	Dorsal margin of the second phalanx	URFXF69	Dorsal margin of the third phalanx	URFXF70	Dorsal margin of the fourth phalanx	URFXF71	Dorsal margin of the fifth phalanx	URFXF72	Dorsal margin of the first phalanx	URFXF73	Dorsal margin of the second phalanx	URFXF74	Dorsal margin of the third phalanx	URFXF75	Dorsal margin of the fourth phalanx	URFXF76	Dorsal margin of the fifth phalanx	URFXF77	Dorsal margin of the first phalanx	URFXF78	Dorsal margin of the second phalanx	URFXF79	Dorsal margin of the third phalanx	URFXF80	Dorsal margin of the fourth phalanx	URFXF81	Dorsal margin of the fifth phalanx	URFXF82	Dorsal margin of the first phalanx	URFXF83	Dorsal margin of the second phalanx	URFXF84	Dorsal margin of the third phalanx	URFXF85	Dorsal margin of the fourth phalanx	URFXF86	Dorsal margin of the fifth phalanx	URFXF87	Dorsal margin of the first phalanx	URFXF88	Dorsal margin of the second phalanx	URFXF89	Dorsal margin of the third phalanx	URFXF90	Dorsal margin of the fourth phalanx	URFXF91	Dorsal margin of the fifth phalanx	URFXF92	Dorsal margin of the first phalanx	URFXF93	Dorsal margin of the second phalanx	URFXF94	Dorsal margin of the third phalanx	URFXF95	Dorsal margin of the fourth phalanx	URFXF96	Dorsal margin of the fifth phalanx	URFXF97	Dorsal margin of the first phalanx	URFXF98	Dorsal margin of the second phalanx	URFXF99	Dorsal margin of the third phalanx	URFXF100	Dorsal margin of the fourth phalanx	URFXF101	Dorsal margin of the fifth phalanx	URFXF102	Dorsal margin of the first phalanx	URFXF103	Dorsal margin of the second phalanx	URFXF104	Dorsal margin of the third phalanx	URFXF105	Dorsal margin of the fourth phalanx	URFXF106	Dorsal margin of the fifth phalanx	URFXF107	Dorsal margin of the first phalanx	URFXF108	Dorsal margin of the second phalanx	URFXF109	Dorsal margin of the third phalanx	URFXF110	Dorsal margin of the fourth phalanx	URFXF111	Dorsal margin of the fifth phalanx	URFXF112	Dorsal margin of the first phalanx	URFXF113	Dorsal margin of the second phalanx	URFXF114	Dorsal margin of the third phalanx	URFXF115	Dorsal margin of the fourth phalanx	URFXF116	Dorsal margin of the fifth phalanx	URFXF117	Dorsal margin of the first phalanx	URFXF118	Dorsal margin of the second phalanx	URFXF119	Dorsal margin of the third phalanx	URFXF120	Dorsal margin of the fourth phalanx	URFXF121	Dorsal margin of the fifth phalanx	URFXF122	Dorsal margin of the first phalanx	URFXF123	Dorsal margin of the second phalanx	URFXF124	Dorsal margin of the third phalanx	URFXF125	Dorsal margin of the fourth phalanx	URFXF126	Dorsal margin of the fifth phalanx	URFXF127	Dorsal margin of the first phalanx	URFXF128	Dorsal margin of the second phalanx	URFXF129	Dorsal margin of the third phalanx	URFXF130	Dorsal margin of the fourth phalanx	URFXF131	Dorsal margin of the fifth phalanx	URFXF132	Dorsal margin of the first phalanx	URFXF133	Dorsal margin of the second phalanx	URFXF134	Dorsal margin of the third phalanx	URFXF135	Dorsal margin of the fourth phalanx	URFXF136	Dorsal margin of the fifth phalanx	URFXF137	Dorsal margin of the first phalanx	URFXF138	Dorsal margin of the second phalanx	URFXF139	Dorsal margin of the third phalanx	URFXF140	Dorsal margin of the fourth phalanx	URFXF141	Dorsal margin of the fifth phalanx	URFXF142	Dorsal margin of the first phalanx	URFXF143	Dorsal margin of the second phalanx	URFXF144	Dorsal margin of the third phalanx	URFXF145	Dorsal margin of the fourth phalanx	URFXF146	Dorsal margin of the fifth phalanx	URFXF147	Dorsal margin of the first phalanx	URFXF148	Dorsal margin of the second phalanx	URFXF149	Dorsal margin of the third phalanx	URFXF150	Dorsal margin of the fourth phalanx	URFXF151	Dorsal margin of the fifth phalanx	URFXF152	Dorsal margin of the first phalanx	URFXF153	Dorsal margin of the second phalanx	URFXF154	Dorsal margin of the third phalanx	URFXF155	Dorsal margin of the fourth phalanx	URFXF156	Dorsal margin of the fifth phalanx	URFXF157	Dorsal margin of the first phalanx	URFXF158	Dorsal margin of the second phalanx	URFXF159	Dorsal margin of the third phalanx	URFXF160	Dorsal margin of the fourth phalanx	URFXF161	Dorsal margin of the fifth phalanx	URFXF162	Dorsal margin of the first phalanx	URFXF163	Dorsal margin of the second phalanx	URFXF164	Dorsal margin of the third phalanx	URFXF165	Dorsal margin of the fourth phalanx	URFXF166	Dorsal margin of the fifth phalanx	URFXF167	Dorsal margin of the first phalanx	URFXF168	Dorsal margin of the second phalanx	URFXF169	Dorsal margin of the third phalanx	URFXF170	Dorsal margin of the fourth phalanx	URFXF171	Dorsal margin of the fifth phalanx	URFXF172	Dorsal margin of the first phalanx	URFXF173	Dorsal margin of the second phalanx	URFXF174	Dorsal margin of the third phalanx	URFXF175	Dorsal margin of the fourth phalanx	URFXF176	Dorsal margin of the fifth phalanx	URFXF177	Dorsal margin of the first phalanx	URFXF178	Dorsal margin of the second phalanx	URFXF179	Dorsal margin of the third phalanx	URFXF180	Dorsal margin of the fourth phalanx	URFXF181	Dorsal margin of the fifth phalanx	URFXF182	Dorsal margin of the first phalanx	URFXF183	Dorsal margin of the second phalanx	URFXF184	Dorsal margin of the third phalanx	URFXF185	Dorsal margin of the fourth phalanx	URFXF186	Dorsal margin of the fifth phalanx	URFXF187	Dorsal margin of the first phalanx	URFXF188	Dorsal margin of the second phalanx	URFXF189	Dorsal margin of the third phalanx	URFXF190	Dorsal margin of the fourth phalanx	URFXF191	Dorsal margin of the fifth phalanx	URFXF192	Dorsal margin of the first phalanx	URFXF193	Dorsal margin of the second phalanx	URFXF194	Dorsal margin of the third phalanx	URFXF195	Dorsal margin of the fourth phalanx	URFXF196	Dorsal margin of the fifth phalanx	URFXF197	Dorsal margin of the first phalanx	URFXF198	Dorsal margin of the second phalanx	URFXF199	Dorsal margin of the third phalanx	URFXF200	Dorsal margin of the fourth phalanx	URFXF201	Dorsal margin of the fifth phalanx	URFXF202	Dorsal margin of the first phalanx	URFXF203	Dorsal margin of the second phalanx	URFXF204	Dorsal margin of the third phalanx	URFXF205	Dorsal margin of the fourth phalanx	URFXF206	Dorsal margin of the fifth phalanx	URFXF207	Dorsal margin of the first phalanx	URFXF208	Dorsal margin of the second phalanx	URFXF209	Dorsal margin of the third phalanx	URFXF210	Dorsal margin of the fourth phalanx	URFXF211	Dorsal margin of the fifth phalanx	URFXF212	Dorsal margin of the first phalanx	URFXF213	Dorsal margin of the second phalanx	URFXF214	Dorsal margin of the third phalanx	URFXF215	Dorsal margin of the fourth phalanx	URFXF216	Dorsal margin of the fifth phalanx	URFXF217	Dorsal margin of the first phalanx	URFXF218	Dorsal margin of the second phalanx	URFXF219	Dorsal margin of the third phalanx	URFXF220	Dorsal margin of the fourth phalanx	URFXF221	Dorsal margin of the fifth phalanx	URFXF222	Dorsal margin of the first phalanx	URFXF223	Dorsal margin of the second phalanx	URFXF224	Dorsal margin of the third phalanx	URFXF225	Dorsal margin of the fourth phalanx	URFXF226	Dorsal margin of the fifth phalanx	URFXF227	Dorsal margin of the first phalanx	URFXF228	Dorsal margin of the second phalanx	URFXF229	Dorsal margin of the third phalanx	URFXF230	Dorsal margin of the fourth phalanx	URFXF231	Dorsal margin of the fifth phalanx	URFXF232	Dorsal margin of the first phalanx	URFXF233	Dorsal margin of the second phalanx	URFXF234	Dorsal margin of the third phalanx	URFXF235	Dorsal margin of the fourth phalanx	URFXF236	Dorsal margin of the fifth phalanx	URFXF237	Dorsal margin of the first phalanx	URFXF238	Dorsal margin of the second phalanx	URFXF239	Dorsal margin of the third phalanx	URFXF240	Dorsal margin of the fourth phalanx	URFXF241	Dorsal margin of the fifth phalanx	URFXF242	Dorsal margin of the first phalanx	URFXF243	Dorsal margin of the second phalanx	URFXF244	Dorsal margin of the third phalanx	URFXF245	Dorsal margin of the fourth phalanx	URFXF246	Dorsal margin of the fifth phalanx	URFXF247	Dorsal margin of the first phalanx	URFXF248	Dorsal margin of the second phalanx	URFXF249	Dorsal margin of the third phalanx	URFXF250	Dorsal margin of the fourth phalanx	URFXF251	Dorsal margin of the fifth phalanx	URFXF252	Dorsal margin of the first phalanx	URFXF253	Dorsal margin of the second phalanx	URFXF254	Dorsal margin of the third phalanx	URFXF255	Dorsal margin of the fourth phalanx	URFXF256	Dorsal margin of the fifth phalanx	URFXF257	Dorsal margin of the first phalanx	URFXF258	Dorsal margin of the second phalanx	URFXF259	Dorsal margin of the third phalanx	URFXF260	Dorsal margin of the fourth phalanx	URFXF261	Dorsal margin of the fifth phalanx	URFXF262	Dorsal margin of the first phalanx	URFXF263	Dorsal margin of the second phalanx	URFXF264	Dorsal margin of the third phalanx	URFXF265	Dorsal margin of the fourth phalanx	URFXF266	Dorsal margin of the fifth phalanx	URFXF267	Dorsal margin of the first phalanx	URFXF268	Dorsal margin of the second phalanx	URFXF269	Dorsal margin of the third phalanx	URFXF270	Dorsal margin of the fourth phalanx	URFXF271	Dorsal margin of the fifth phalanx	URFXF272	Dorsal margin of the first phalanx	URFXF273	Dorsal margin of the second phalanx	URFXF274	Dorsal margin of the third phalanx	URFXF275	Dorsal margin of the fourth phalanx	URFXF276	Dorsal margin of the fifth phalanx	URFXF277	Dorsal margin of the first phalanx	URFXF278	Dorsal margin of the second phalanx	URFXF279	Dorsal margin of the third phalanx	URFXF280	Dorsal margin of the fourth phalanx	URFXF281	Dorsal margin of the fifth phalanx	URFXF282	Dorsal margin of the first phalanx	URFXF283	Dorsal margin of the second phalanx	URFXF284	Dorsal margin of the third phalanx	URFXF285	Dorsal margin of the fourth phalanx	URFXF286	Dorsal margin of the fifth phalanx	URFXF287	Dorsal margin of the first phalanx	URFXF288	Dorsal margin of the second phalanx	URFXF289	Dorsal margin of the third phalanx	URFXF290	Dorsal margin of the fourth phalanx	URFXF291	Dorsal margin of the fifth phalanx	URFXF292	Dorsal margin of the first phalanx	URFXF293	Dorsal margin of the second phalanx	URFXF294	Dorsal margin of the third phalanx	URFXF295	Dorsal margin of the fourth phalanx	URFXF296	Dorsal margin of the fifth phalanx	URFXF297	Dorsal margin of the first phalanx	URFXF298	Dorsal margin of the second phalanx	URFXF299	Dorsal margin of the third phalanx	URFXF300	Dorsal margin of the fourth phalanx	URFXF301	Dorsal margin of the fifth phalanx	URFXF302	Dorsal margin of the first phalanx	URFXF303	Dorsal margin of the second phalanx	URFXF304	Dorsal margin of the third phalanx	URFXF305	Dorsal margin of the fourth phalanx	URFXF306	Dorsal margin of the fifth phalanx	URFXF307	Dorsal margin of the first phalanx	URFXF308	Dorsal margin of the second phalanx	URFXF309	Dorsal margin of the third phalanx	URFXF310	Dorsal margin of the fourth phalanx	URFXF311	Dorsal margin of the fifth phalanx	URFXF312	Dorsal margin of the first phalanx	URFXF313	Dorsal margin of the second phalanx	URFXF314	Dorsal margin of the third phalanx	URFXF315	Dorsal margin of the fourth phalanx	URFXF316	Dorsal margin of the fifth phalanx	URFXF317	Dorsal margin of the first phalanx	URFXF318	Dorsal margin of the second phalanx	URFXF319	Dorsal margin of the third phalanx	URFXF320	Dorsal margin of the fourth phalanx	URFXF321	Dorsal margin of the fifth phalanx	URFXF322	Dorsal margin of the first phalanx	URFXF323	Dorsal margin of the second phalanx	URFXF324	Dorsal margin of the third phalanx	URFXF325	Dorsal margin of the fourth phalanx	URFXF326	Dorsal margin of the fifth phalanx	URFXF327	Dorsal margin of the first phalanx	URFXF328	Dorsal margin of the second phalanx	URFXF329	Dorsal margin of the third phalanx	URFXF330	Dorsal margin of the fourth phalanx	URFXF331	Dorsal margin of the fifth phalanx	URFXF332	Dorsal margin of the first phalanx	URFXF333	Dorsal margin of the second phalanx	URFXF334	Dorsal margin of the third phalanx	URFXF335	Dorsal margin of the fourth phalanx	URFXF336	Dorsal margin of the fifth phalanx	URFXF337	Dorsal margin of the first phalanx	URFXF338	Dorsal margin of the second phalanx	URFXF339	Dorsal margin of the third phalanx	URFXF340	Dorsal margin of the fourth phalanx	URFXF341	Dorsal margin of the fifth phalanx	URFXF342	Dorsal margin of the first phalanx	URFXF343	Dorsal margin of the second phalanx	URFXF344	Dorsal margin of the third phalanx	URFXF345	Dorsal margin of the fourth phalanx	URFXF346	Dorsal margin of the fifth phalanx	URFXF347	Dorsal margin of the first phalanx	URFXF348	Dorsal margin of the second phalanx	URFXF349	Dorsal margin of the third phalanx	URFXF350	Dorsal margin of the fourth phalanx	URFXF351	Dorsal margin of the fifth phalanx	URFXF352	Dorsal margin of the first phalanx	URFXF353	Dorsal margin of the second phalanx	URFXF354	Dorsal margin of the third phalanx	URFXF355	Dorsal margin of the fourth phalanx	URFXF356	Dorsal margin of the fifth phalanx	URFXF357	Dorsal margin of the first phalanx	URFXF358	Dorsal margin of the second phalanx	URFXF359	Dorsal margin of the third phalanx	URFXF360	Dorsal margin of the fourth phalanx	URFXF361	Dorsal margin of the fifth phalanx	URFXF362	Dorsal margin of the first phalanx	URFXF363	Dorsal margin of the second phalanx	URFXF364	Dorsal margin of the third phalanx	URFXF365	Dorsal margin of the fourth phalanx	URFXF366	Dorsal margin of the fifth phalanx	URFXF367	Dorsal margin of the first phalanx	URFXF368	Dorsal margin of the second phalanx	URFXF369	Dorsal margin of the third phalanx	URFXF370	Dorsal margin of the fourth phalanx	URFXF371	Dorsal margin of the fifth phalanx	URFXF372	Dorsal margin of the first phalanx	URFXF373	Dorsal margin of the second phalanx	URFXF374	Dorsal margin of the third phalanx	URFXF375	Dorsal margin of the fourth phalanx	URFXF376	Dorsal margin of the fifth phalanx	URFXF377	Dorsal margin of the first phalanx	URFXF378	Dorsal margin of the second phalanx	URFXF379	Dorsal margin of the third phalanx	URFXF380	Dorsal margin of the fourth phalanx	URFXF381	Dorsal margin of the fifth phalanx	URFXF382	Dorsal margin of the first phalanx	URFXF383	Dorsal margin of the second phalanx	URFXF384	Dorsal margin of the third phalanx	URFXF385	Dorsal margin of the fourth phalanx	URFXF386	Dorsal margin of the fifth phalanx	URFXF387	Dorsal margin of the first phalanx	URFXF388	Dorsal margin of the second phalanx	URFXF389	Dorsal margin of the third phalanx	URFXF390	Dorsal margin of the fourth phalanx	URFXF391	Dorsal margin of the fifth phalanx	URFXF392	Dorsal margin of the first phalanx	URFXF393	Dorsal margin of the second phalanx	URFXF394	Dorsal margin of the third phalanx	URFXF395	Dorsal margin of the fourth phalanx	URFXF396	Dorsal margin of the fifth phalanx	URFX
--	----------------	-------------	-------------------	--------	---	--------	--	--------	--	--------	---	--------	---	--------	--------------------------------------	--------	--	--------	-------------------------	--------	---------------------------------	---------	-----------------------------	---------	------------------------------	---------	--	---------	---	---------	--	---------	---	---------	--	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	---------	--	---------	---	---------	---	---------	--	---------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	----------	---	----------	--	----------	---	----------	--	----------	---	------

References

1. ↑ [1.00 1.01 1.02 1.03 1.04 1.05 1.06 1.07 1.08 1.09 1.10 1.11 1.12 1.13 1.14 1.15 1.16 1.17 1.18 1.19 1.20 1.21 1.22 1.23 1.24 1.25 1.26 1.27 1.28 1.29 1.30 1.31 1.32 1.33 1.34](#) Serge van Sint Jan “[Color Atlas of Skeletal Landmark Definitions: Guidelines for Reproducible Manual and Virtual Palpations](#)” 2007 - Churchill Livingstone
2. ↑ [2.0 2.1](#) Leardini, A., M.G. Benedetti, L. Berti, D. Bettinelli, R. Natio, and S. Giannini. “[Rear-foot, Mid-foot and Fore-foot Motion during the Stance Phase of Gait.](#)” *Gait & Posture* 25 (2007): 453-55

From: <https://has-motion.com/wiki/> - **Wiki Documentation Site**

Permanent link:
https://has-motion.com/wiki/doku.php?id=other:ior_gait:documentation:marker_sets&rev=1718717127

Last update: **2024/06/18 13:25**

