

RT_Plugin_MAC_Cortex

Overview

The MAC Cortex™ Plugin is used in [Visual3D Server™](#), and Visual3D RT™, it connects to [Motion Analysis™](#) software.

Plugin Interface



There are four sets of tabs on the left hand side of the screen. These tabs are [Hardware](#), [Signals](#), [C3D Params](#), and [Force Plates](#). Each of these tabs controls a different portion of the plugins functionality. On the right hand side of the window is a [3D view screen](#) which shows a graphical representation of the data the server is receiving from the MAC Cortex hardware.



The “Plugin Help” button leads to this web page should you need to reference this page in the future.



The “About Plugin” button has details on the plugin being used including version, release date, copy right, support contact information and the hardware SDK version.

Frame Info



Current frame The current frame number on the top left hand side on the screen (where it is currently blank) keeps a count of the current frame which is assigned by the hardware or software system.

Missed frames The missed frames number is on the bottom left (where it is currently blank). Whenever there is a gap in frames sent by the hardware or software system a blank frame is inserted. The dropped frames count keeps a record of both the number of frames which are blank.

Hardware time The hardware time number is on the top, in the centre (where it is currently blank). This time stamp is created by the hardware being used and is how long, in milliseconds, the hardware has been running.

Server time The server time number is on the bottom, in the centre (where it is currently blank). This time stamp is created by the server and is how long, in milliseconds, it has been connected to the hardware or software which is streaming to it.

Hardware latency The hardware latency is on the top right hand side (where it is currently blank). This indicated the delay created by the hardware system.

Server latency The hardware latency is on the top right hand side (where it is currently blank). This indicates any delay introduced by the server.

Hardware

== IP Address ==

In order to connect with the data that is being broadcast for the server the ServerIP address must be entered into the dialogue box at the top of the plugin tab. The IP address used in the image above is 192.168.1.102. MAC Cortex™ does not allow for the “localhost” or 127.0.0.1 shortcuts to be used. You will need to use your actual IP address. See [Trouble Shooting](#) for how to check your IP address.

== Calibration File ==

The Cal file box is for systems which use force plates. The calibration file indicates where the force platforms are located relative to the lab coordinate system. This file is required in order for force data to be included in the graphics and in the data which is streamed to Visual3D server. Consult your MAC Cortex™ documentation for instructions on how to generate a calibration file. You should generate and keep a calibration file every time that the force plates may have been moved. Files can be selected using the Browse button and can be removed using the Remove button.

== Gen Scale ==

The bits and volts drop down menus tell the server how to process the force data that it receives. If you insert the wrong bits or volts values your calculations and data will be erroneous. The bits and volts were determined when the force platforms were set up; see your documentation or the person responsible for the set up to get these values.

Signals

✖

The table in the signals tab holds all of the numerical data from various sources which is being streamed to the program. This data may include, but is not limited to, marker co-ordinates, force data, EMG readings, accelerometer data, etc. The first column entitled “Label” has the number or name which has been assigned to the source of the data. The following four columns are all entitled “Value”; they have the numerical data associated with the source. For markers it is the x, y, and z, coordinates and then the residual. For force data there is only one value associated, which is the magnitude of the force.

✖

The signals graph allows you to see a graphical representation of the values in the table. Selecting one of the rows in the table will cause the graph to record the values in real time. The x axis has the frame numbers and the y axis has the numbers in the value column. If more than one column has data there will be multiple lines drawn. The graph uses the MKS system of units.

C3D Params

✖

When this tab is viewed in Visual3D Server it gives additional details regarding the data that is being collected by the server and streamed to a client. This allows you to check details such as force plate types and locations.

✖

The top table contains a list of the parameters and their subsections. By expanding the parameters and selecting one of the subsections the data from the subsection will be presented in the table at the bottom of the tab.

Force Platform

The minimum force management box is for screening out unwanted noise from the force platforms. The default value is 20.00 N meaning that any force equal to or lower than 50.00 N will not be displayed by the system. It does not affect forces that are greater than the specified limit. The minimum force may be set at any value between 0 and 9,999.90 N.

The baseline force measurement tool allows you to determine the average amount of noise present in your force platform system and remove it from the forces displayed by the system and sent to any client programs such as a biofeedback software. The baseline frame number is set at 10 which means that the baseline function will take a sample of 10 consecutive frames. The frame number may be set at any whole number from 1 to 9,999. When you click the "Measure Baseline" button it will take a sample of frames (as many or as little as indicated by the baseline frame number box.) The average force will be calculated for these frames. This value will subsequently be subtracted from all of the following frames. The graphics displayed and the data passed on to any client programs will be the initial value minus the calculated average. When the "Zero Baseline" button is pressed the forces will no longer have the baseline number subtracted and the forces displayed and given to any client programs will be the value taken from the force plate.

Both force management tools may be used at the same time.

The "Use force structure" checkbox controls the force structure, which is the amalgamation of multiple forces. If the checkbox is not selected then the force structure will not be calculated. This means that the force structure will not appear in the graphics and that the data for a force structure will not be available for any client programs.

3D View

The right hand side of the screen provides a real-time 3D graphical representation of the data being streamed.

Save 3D View



This button allows you to save what is currently being displayed in the 3D view graphics area. The image can be saved in a number of standard formats, such as .bmp, .ico, .jpeg, .jpg, .pbm, .pgm, .png, .ppm, .tif, .tiff, .wbmp, .xbm, or .xpm. This function is ideal for producing graphics for papers or presentations. Saved images can be edited with 3rd party graphics programs if needed.

Copy 3D View to clipboard



This button allows you to copy what is currently being displayed in the 3D view graphics area to the clipboard. This allows you to paste the image into any program that supports copy and paste. This is useful for working with presentation or graphics editing software.

Reset camera 3D View



This button brings the 3D viewer back to it's beginning position. You may find that having used the rotation, translation, and zooming capabilities of the 3D view you have lost track of where you are in the graphics, which can make returning to a proper view of your data challenging. The reset view button removes this frustration and reorients you.

3D View Preferences



This button allows you to change how the graphics that are presented to you. Changes happen in real time. At the bottom is the “Restore Defaults” button which resets all of the options to their standard setting.

= Scene =



Lighting on - When selected graphics that are suppose to be 3 dimensional have shadows to help simulate shape. When not selected objects are one solid color.

Reset camera view on connect - When this box is selected and you connect or reconnect to hardware the camera view will reset to it's default.

Alpha-blending on - When this box is selected the graphics are slightly transparent and allow you to see what is behind them

Anti-ailiasing on - When this box is selected all the lines in the graphics are smooth. If it is not selected the lines that are not either horizontal or vertical change directions portion by portion.

= Background =

Color - This allows you to change the background color to anyone you chose.

= Floor =

When the check box next to the floor is not checked the floor will not appear in the 3D View. In this case all the bellow options will be greyed out as they will have no effect if the floor is not present.

X-size - This allows you to customize how many panels the floor is composed of in the x direction. the minimum is 1 and maximum is 100. The default size is 5.00.

Y-size - This allows you to customize how many panels the floor is composed of in the y direction. the minimum is 1 and maximum is 100. The default size is 5.00.

X-offset - This allows you to shift the entire floor along the x-axis either in the positive or negative direction. The maximum value allowed is 100 and the minimum value is -100. The default setting is 3.00.

Y-offset - This allows you to shift the entire floor along the y-axis either in the positive or negative direction. The maximum value allowed is 100 and the minimum value is -100. The default setting is 3.00.

Z-offset - This allows you to shift the entire floor along the z-axis either in the positive or negative

direction. The maximum value allowed is 100 and the minimum value is -100. The default setting is -0.01.

Color - This allows you to change the floor color to anyone you chose.

Show grid - This allows you to remove the grid lines from the floor by unchecking the checkbox.

Grid line width - This allows you to change the thickness of the gridlines. The minimum value is 1 and the maximum value is 10. The default width is 2.00.

Grid color - This allows you to change the grid line color to anyone you chose.

= Axes =

When the check box next to the axis is not checked the axis will not appear in the 3D View. In this case all the bellow options will be greyed out as they will have no effect if the axis is not present.

Line width - This allows you to change the thickness of the axis lines. The minimum value is 1 and the maximum value is 10. The default value is 2.00.

X-Axis Color - This allows you to change the x-axis line color to anyone you chose.

Y-Axis Color - This allows you to change the y-axis line color to anyone you chose.

Z-Axis Color - This allows you to change the z-axis line color to anyone you chose.

= Force Plates =

When the check box next to the force plates is not checked the force plates will not appear in the 3D View. In this case all the bellow options will be greyed out as they will have no effect if the force plates are not present.

Plate Color - This allows you to change the force plate color to anyone you chose.

Show labels - This allows you to remove the force plate label, (such as FP1, FP2 . . . etc) from the 3D view.

Label Color - This allows you to change the force plate label color to anyone you chose.

Show forces - When checked the forces will be show in the 3D view using arrows.

Force Type A Color - This allows you to change the color of the raw force arrows to anyone you chose.

Force Type B Color - This allows you to change the color of the force structure arrows to anyone you chose.

= Tracking Markers =

When the check box next to the tracking markers is not checked the tracking markers will not appear in the 3D View. In this case all the bellow options will be greyed out as they will have no effect if the tracking markers are not present.

Radius - This allows you to change the radius of the markers. The maximum value is 0.100 the minimum value is 0.001. The default radius is 0.020.

Marker Color - This allows you to change the color of the marker color to anyone you chose.

Show Labels - This allows you to remove the marker labels from the 3D view.

Label Color - This allows you to change the color of the label to anyone you chose.

= Segment Markers =



When the check box next to the segment markers is not checked the segment markers will not appear in the 3D View. In this case all the bellow options will be greyed out as they will have no effect if the segment markers are not present.

Radius - This allows you to change the radius of the markers. The maximum value is 0.100 the minimum value is 0.001. The default radius is 0.020.

Marker Color - This allows you to change the color of the marker color to anyone you chose.

Show Labels - This allows you to remove the marker labels from the 3D view.

Label Color - This allows you to change the color of the label to anyone you chose.

= Segment Lines =

When the check box next to the segment lines is not checked the segment lines will not appear in the 3D View. In this case all the bellow options will be greyed out as they will have no effect if the segment lines are not present.

Line Width - This allows you to change the thickness of the segment lines. The minimum value is 1 and the maximum value is 10. The default value is 1.00.

Color - This allows you to change the color of the segment lines to anyone you chose.

= Segment Geometry =

When the check box next to the segment geometry is not checked the segment geometry will not appear in the 3D View. In this case all the bellow options will be greyed out as they will have no effect if the segment geometry is not present.

Color - This allows you to change the color of the segment geometry to anyone you chose.

= Segment Bones =

When the check box next to the segment bones is not checked the segment bones will not appear in the 3D View. In this case all the bellow options will be greyed out as they will have no effect if the segment bones are not present.

Color - This allows you to change the color of the segment bones to anyone you chose.

From:
<https://has-motion.com/wiki/> - Wiki Documentation Site

Permanent link:
https://has-motion.com/wiki/doku.php?id=other:visual3dserver:plugins:rt_plugin_mac_cortex&rev=1721157796

Last update: **2024/07/16 19:23**

