

Introduction

How to Modify Explicit Settings for Report Graphs

By default, the Y axis of a graph will be graphed from max to min:

[ReportGraphExplicit_2.gif](#){ [{/images/6/67/ReportGraphExplicit_2.gif](#) **Changing the settings will allow you to manually specify the max/min value:**

[ReportGraphExplicit_5.gif](#)}}

Consider a graph created with the settings on the right:

Since the Y-Axis Scale is set to Global Min to Max, the graph that is created has the range of data for the y-axis determined by the maximum and minimum values of the signals graphed:

[ReportGraphExplicit_2.gif](#){ [{/images/6/67/ReportGraphExplicit_2.gif](#) [ReportGraphExplicit_1.png](#)}}

To change the range:

1. In the Reports tab, select the graph using the Left Mouse Button - Handles will appear indicating that the graph is active - If the handles don't appear, the graph is not selected!! 2. Right click on the selected graph 3. Select Format Graph	ReportGraphExplicit_3.gif
--	--

4. Set the y-axis range: - Check Modify Y Axis Scale - Specify the Y Min/Y Max 5. Click Apply	ReportGraphExplicit_4.png
---	--

The graph has now been explicitly set to these properties: [ReportGraphExplicit_5.gif](#)

The Report Graph dialog will now reflect that the settings are explicit. If the graph template is saved, these explicit settings will be saved in the Report Template	ReportGraphExplicit_6.png
---	--

From:
<https://has-motion.com/wiki/> - **Wiki Documentation Site**

Permanent link:
https://has-motion.com/wiki/doku.php?id=visual3d:documentation:reports:explicit_settings_for_report_graphs&rev=1718648613

Last update: **2024/06/17 18:23**

